## BIOMARK Laboratories-INDIA www.biomarklabs.com TECHNICAL SHEET

BF055	FRASER ENRICHMENT SUPPLEMENT	
FORMULA		
Ingredients : (Per vial, sufficient for 500 ml r	iedium)	
Acriflavine hydrochloride Nalidixic acid	25mg 4mg	
Directions :		
water. Mix well and a	s of one vial aseptically with 10 ml of sterile distilled eptically add it to 990 ml sterile, cooled (45-50°C) chment Broth Base (B1045). Mix well and dispense a	ЭS
Precautions : 1. For Laboratory Use 2. Follow proper, esta disposing of infection	plished laboratory procedures in handling and	
	nt recommended for enrichment of Listeria s and environmental specimens.	
Storage:Between 2 to 8°C.		
Packing :5 vial		